



CORA STYLES LAB SUPPLIES, LLC

PO Box 431
TALENT OR 97540
USA

PHONE (541) 535-6775
FAX (678) 693-5704

YEAST REPLICATOR BLOCKS WITH COLLARS

Both the 4-inch and the 6-inch aluminum replicating blocks are machined to have the same space between the inside ring of the collar and the lower ring of the replicating block. When a velveteen square is placed on the block and the collar pressed down to hold it in place, it may seem that the 6-inch block doesn't grip the cloth as tightly as the 4-inch block does. This is because the cloth bunches up more on the block with the smaller diameter.

If you wish the velveteen to be more tightly gripped on either of the blocks, we recommend wrapping 1/2- or 1/4-inch wide masking tape around the lower ring once or more.